

Available Configuration

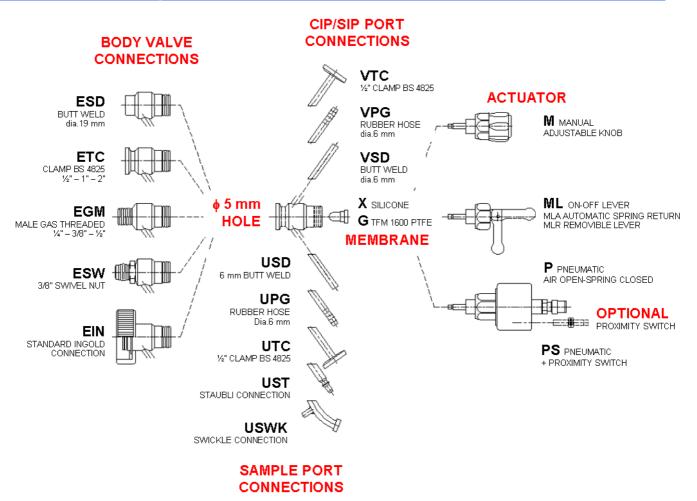

Ordering information

To specify the part completely, start with the product description and select the additional options as shown below :

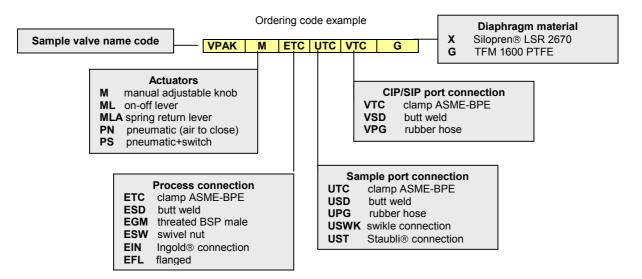

VPAK - PTFE Diaphragm

Technical specification

TFM 1600 PTFE diaphragm are obtained, by means of machining on CCN, from **TFM™ 1600 PTFE** bar, Poly-Tetra-Fluor-Ethylene from Dyneon[™]- 3M. compound are manufactured conforming to the following norms :

- FDA Code of Federal Regulation
- CFR under title 21, paragraph 177.1550
- U.S.Pharmacopoeia Class VI

Maintenance :

TFM 1600 PTFE diaphragm must be replaced once every year, but in the event of intensive use, sterilisation and cleaning it may be necessary to replace it more frequently.


AR-Inox guaranties one year of lasting by normal

use means (silicone has to be changed approx every 4-5 months)

Temperature......121°C (150°C for short period only) Steam pressure......1,5 up to 2 bar max.

Sample.....1-5 a day

Chemical resistance : Is not attacked by common chemicals

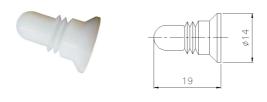

TFM 1600 PTFE				
Ordering code		MVPAK-G		
Material		TFM 1600 PTFE		
Colour		White		
Powder compound properties				
Property	Value	Unit	Test Method	
Bulk density	830	g/l	ASTM D 4894-98a	
Average particle size	450	μ	ASTM D 4894-98a	
Mechanical properties, measured at 23°C (73°F)				
Tensile Strength	4600	psi	ASTM D 4894-98a	
Elongation at break	450	%	ASTM D 4894-98a	
Specific gravity	2.16	g/cc	ASTM D 4894-98a	
Shrinkage	3.5	%	ASTM D 4894-98a	
Tensile Modulus	94,250	psi	ASTM D 638	
Deformation under Load		%	ASTM D 621	
2175 psi – 24 hrs	8			
2175 psi – 100 hrs	9			
2175 psi – permanent	4			
Thermal properties				
Flammability	V-0		UL94	
Melt point (initial)	342 ±10	°C	ASTM D 4894-98a	

VPAK - SILICONE Diaphragm

Technical specification

Build by **Silopren®** LSR 2670 a silicone elastomer composed from two components, it is supplied by GE Bayer Silicones and it is obtained with a modern system of injection molding.

After molding, they under go a heat treatment (post-cured) of 4 hours at 200°C inside a furnace with forced ventilation.

All diaphragms are manufactured in 100% silicone of medical grade conforming to the following norms:

- FDA Code of Federal Regulation
- CFR 177.2600
- U.S. Pharmacopeia Class VI


Maintenance :

Average live of a silicone diaphragm is 4-5 months of lasting by normal use means

Temperature max......121°C Steam pressure......1,5 up to 2 bar max. Process pressure.......6 bar max. Sample......1-5 a day

not recommended for use with alcool, solvent, glicerina

Silopren LSR 2670			
Ordering code	MVPAK-X		
Material	Silopren® LSR 2670		
Colour	White		
Hardness	60 Shore		
Tensile strength	10,5 MPa		
Wear resistance	Less suitable		
Tear resistance	Very good		
Resistance to Weather and Ozone	Excellent		
Resistance to Hydrolysis (water and steam)	Good		
Resistance to Chemicals (acids / bases)	Suitable		
Resistance to mineral oil and gas	Less suitable		
Air and Gas density	Not suitable		

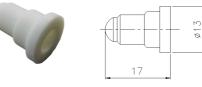

WARNING !

THE INSPECTION CHECK AND MAINTENANCE INTERVALS CAN DEPENDS FROM OPERATING CONDITIONS AS TEMPERATURE, PRESSURE AND FLUID TYPE HOWEVER, ALWAYS REMAIN TO THE USER, THE RESPONSABILITY TO FIX A SUITABLE PROCEDURE WITH THE INTERVALS AND FORMALITY IN ORDER TO REPLACE THE DIAPHRAGM BASED ON OWN EXPERIENCE.